ISENO3181

Basic User Manual

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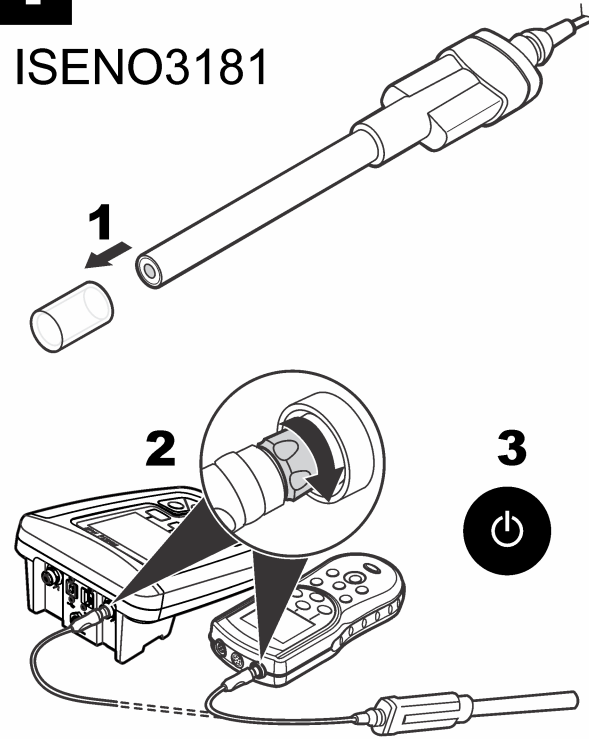
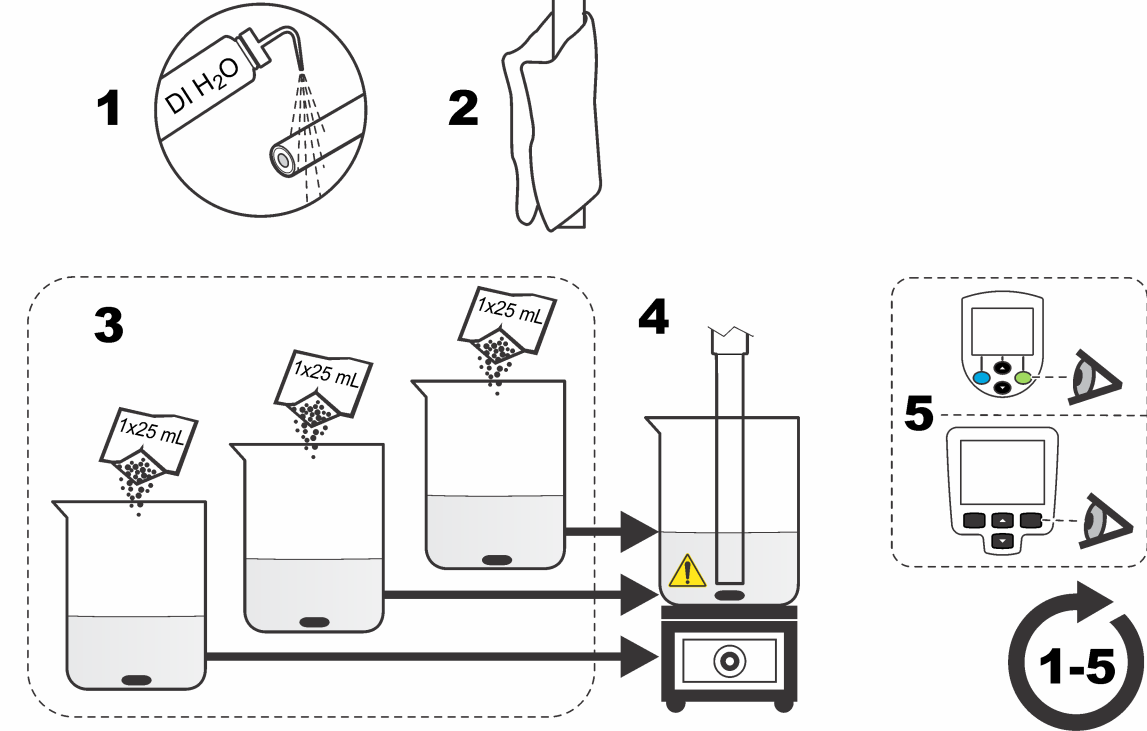
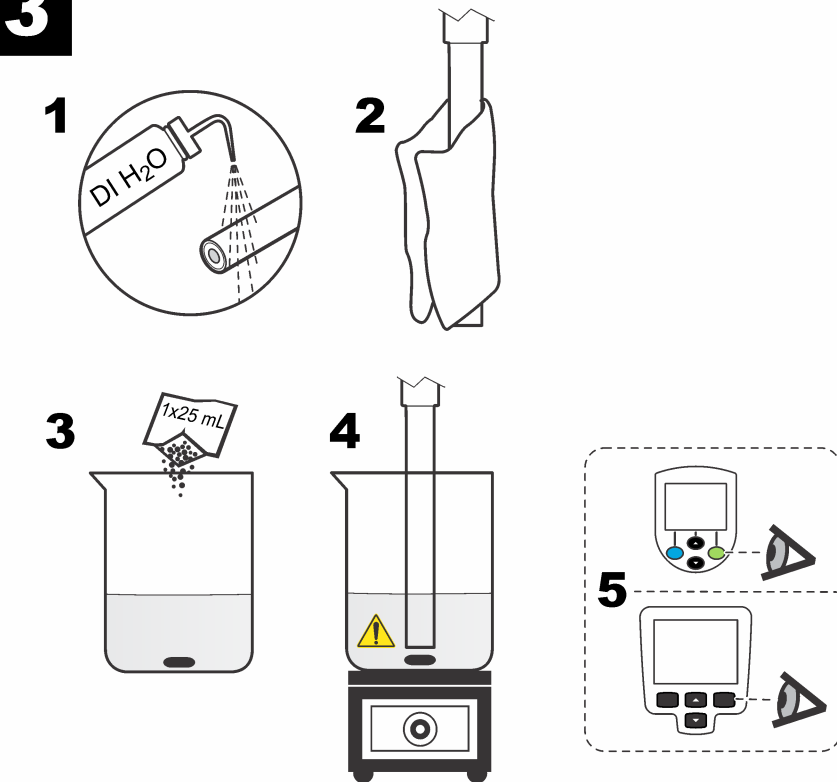
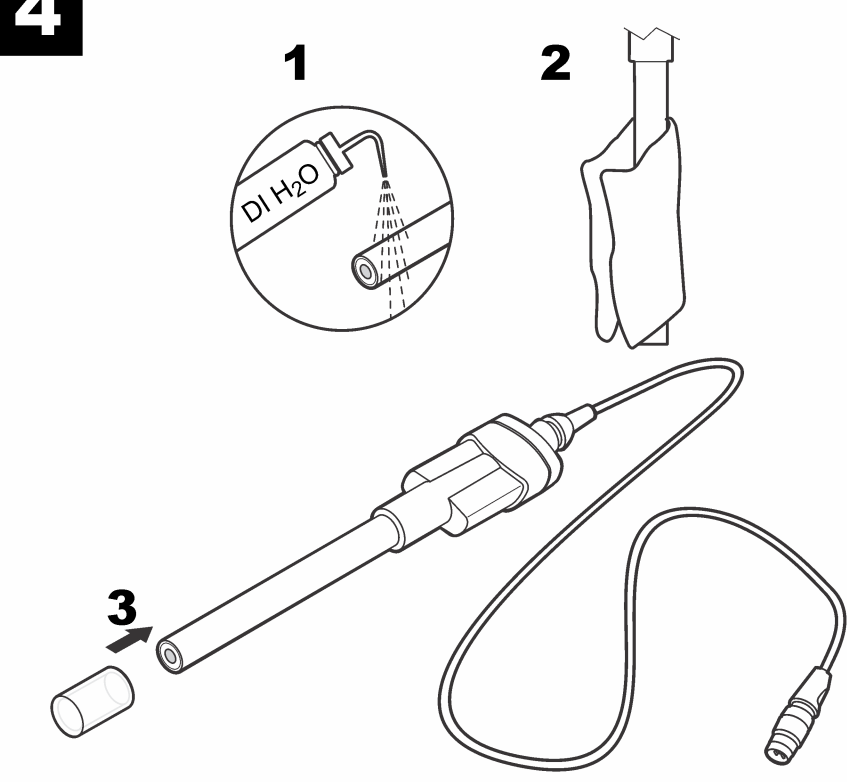
www.hach.com

1**ISENO3181****2****3****4****EN****ISENO3181 probe**

Additional information is available on the manufacturer's website.

Introduction

The probe is a digital combination probe with a non-refillable double junction reference and a built-in temperature sensor.

SAFETY NOTE

Before use with hazardous substances, find out the relevant hazards, protective measures and what to do in an emergency.

Collect the following items:

Nitrate standard solutions and ionic strength adjusters; beakers, 50 mL; stirrer plate; DI water; stir bars; lint-free cloth.

1 Connect the probe**2 Calibration**

▲ Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.

3 Measurement (direct method)

Temperatures of calibration standards and samples should be kept within ± 2 °C of each other for optimal results.

4 Storage**Technical information**

Range	0.1 mg/L (7×10^{-6} M) to 14,000 mg/L (1 M) nitrate nitrogen
pH Range	pH 2 to 11, must be adjusted to 3 to 5 by nitrate ISA
Operating temperature	5 to 50 °C (41 to 122 °F)
Minimum sample volume	25 mL
Ionic strength adjusters	Hach nitrate ISA (1 powder pillow per 25 mL standard or sample)
Default standard solutions	1, 10 and 100 mg/L Hach nitrate nitrogen standard solutions

Maintenance and storage

The probe can be stored dry. To protect the sensing element, rinse with DI water and blot dry with a lint-free cloth. Install the sensor protection cap.

Warranty

1 year for manufacturing faults. This warranty does not cover improper use or wear.

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