50 52 T

User Manual

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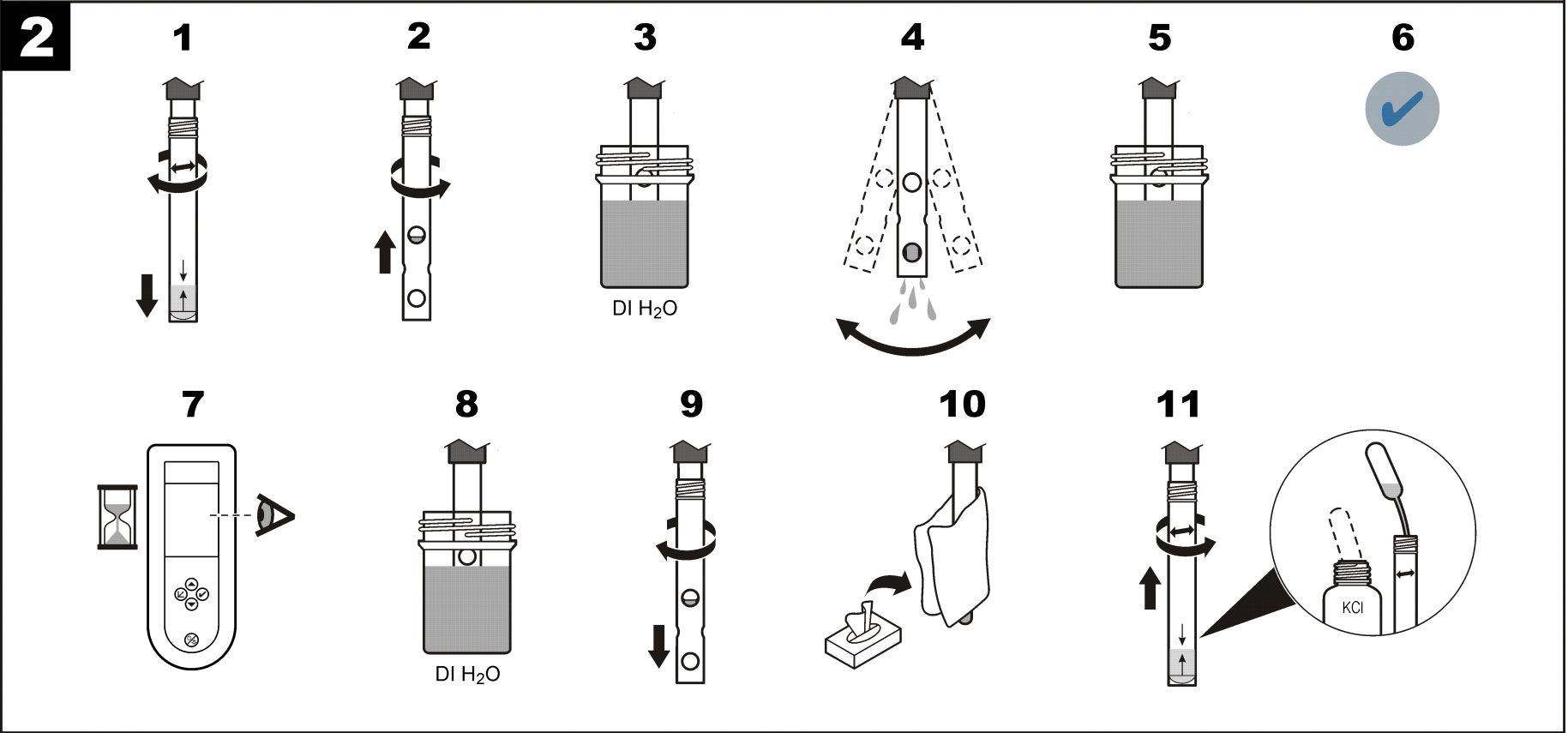
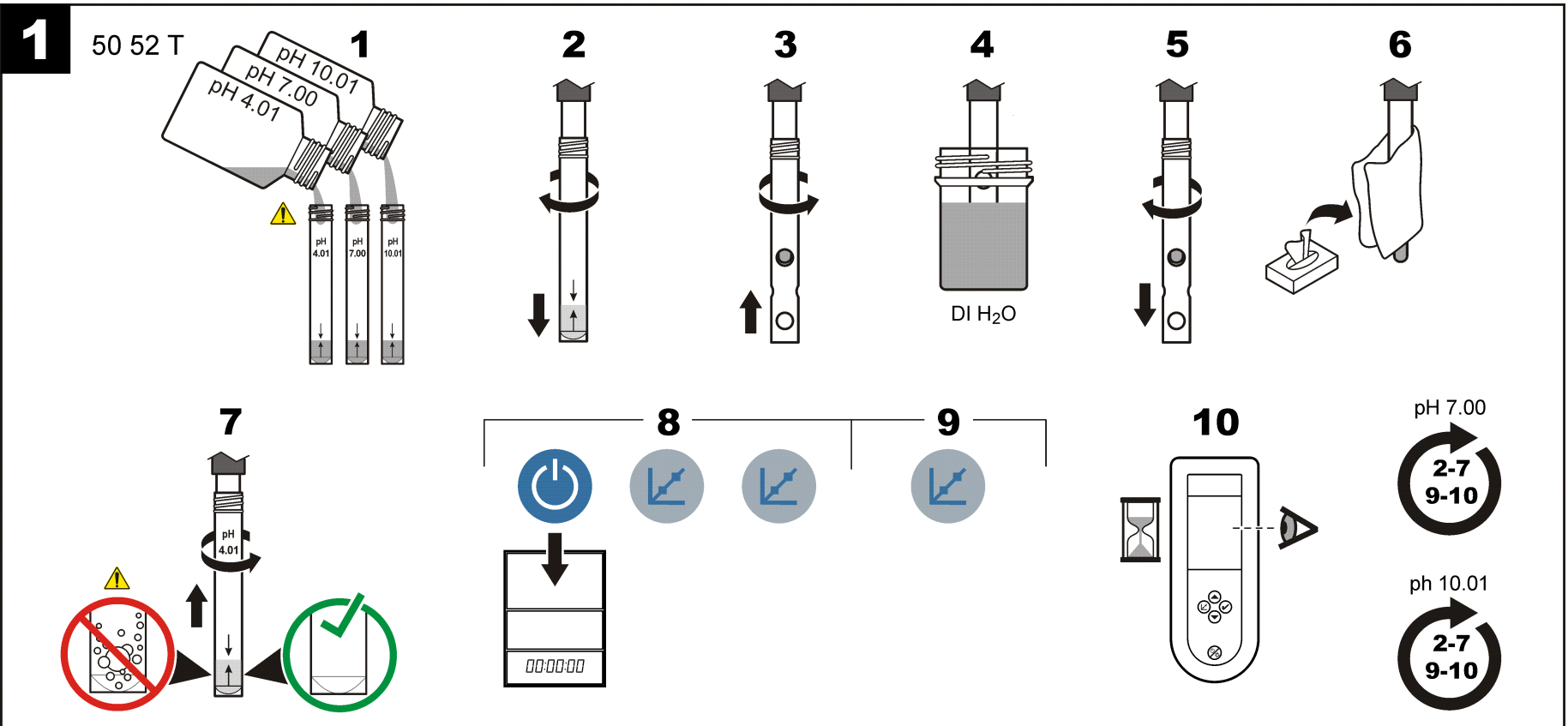
EN

50 52 T probe
Refer to the meter manual for more information.

Caution
 ⚠ *Personal injury hazard. Glass components can break. Handle with care to prevent cuts.*
 ⚠ *Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.*

Introduction
This probe is a combination gel-filled electrode with a built-in temperature sensor for pH measurements in low-conductivity, dirty, viscous and high temperature samples.

Technical information

Range	0 to 14 pH
Junction	PTFE
Reference	Ag/AgCl
Temperature	Pt 1000
Minimum depth	12 mm (0.5 in.)
Operating temp	0–100 °C (32–212 °F)
Connector type	MP-5

Preparation
Install the protective shroud.

Connect the probe
Caution
⚠ *Do not turn or twist the probe connector.*

1 Calibration
Caution
⚠ *Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.*
⚠ *Rinse sensor thoroughly between standards.*

For 1 or 2-point pH calibration: push ✓ after 1 or 2 standards.

2 Sample measurement
Continuous: push ✓ 2 times.
Change the parameter: push and hold ✓ during stabilization.

Maintenance and storage
Cleaning: soak the probe in a cleaning solution. Rinse with deionized water.
Hydration: soak the dry probe in a dilute HCl solution for several hours.
Storage: put the probe in the protector with KCl 3 M solution.

Troubleshooting
Be sure that the correct probe type is used. Do not hold the tubes by hand to prevent temperature changes.

Warranty
6 months for manufacturing faults only. Damage from use is not covered.

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