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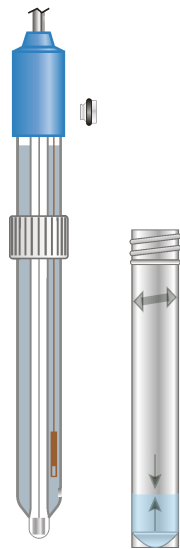
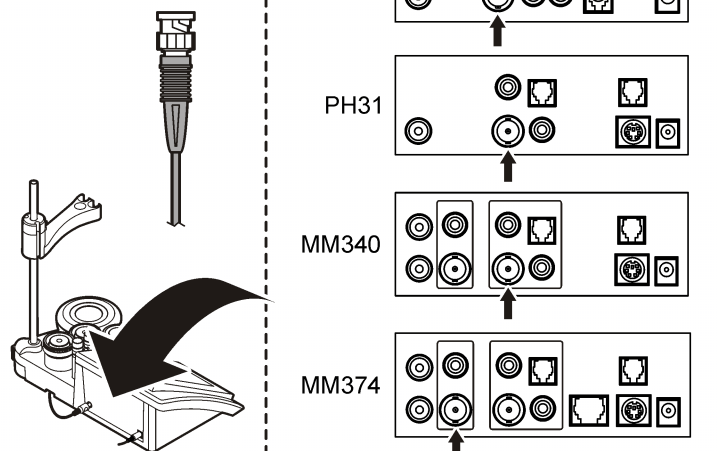
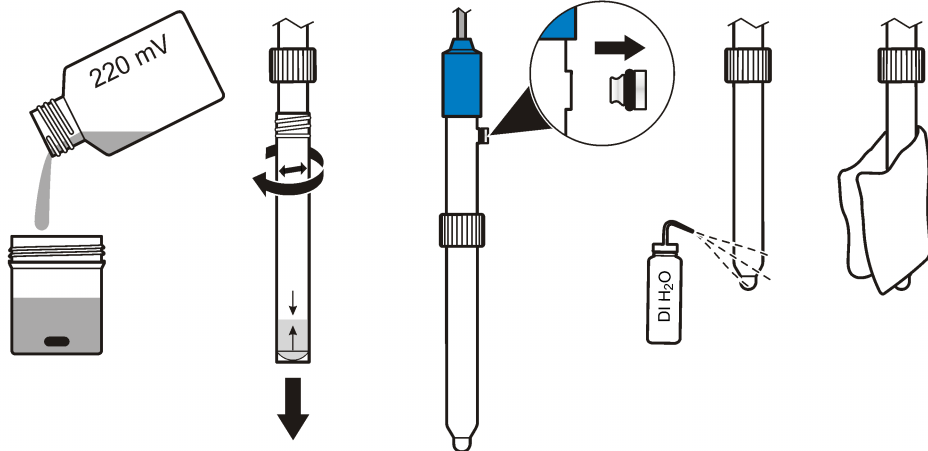
1 ⚠

50 57

Refer to the meter manual for more information.

**Caution**

⚠ *Personal injury hazard. Glass components can break. Handle with care to prevent cuts.*

⚠ *Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.*

**Introduction**

This refillable probe is a platinum electrode with a plate sensor for oxidation reduction potential (ORP) measurement.

**Technical information**

Range	± 2000 mV
Indicator	Platinum
Junction	Ceramic
Reference	Ag/AgCl
Minimum depth	12 mm (0.5 in.)
Operating temp	0 to 80 °C (32 to 176 °F)
Limits	Colloids or solids in suspension; viscous or dirty solutions

**Preparation**

Remove the storage protector from the probe.

**1 Connect the probe****2 Calibration**

⚠ *Air bubbles in or under the probe tip when submerged can cause slow stabilization or error in measurement.*

**3 Measurement****Maintenance and storage**

Cleaning: rinse the junction with distilled water after measurement. Use a mild abrasive such as toothpaste to clean the platinum sensor.

Refill the electrode with electrolyte as needed.

Storage: put the probe in the protector with KCl 3 M solution. Be sure that the refill port is closed.

**Troubleshooting**

Be sure that the samples are suitable for the application.

Keep the junction and platinum sensor free of dirt.

**Warranty**

1 year for manufacturing faults only. Damage from use is not covered.

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