



INSTRUCTION SHEET

Odyssey DR/2500 Instrument Update Module

Using an update module, the Odyssey DR/2500 Spectrophotometer can be updated in the field to include the latest versions of chemistry methods or software from Hach Company. This instruction sheet covers the use of two types of Odyssey Instrument Update Modules.

1. The Hach Program Update Module, Cat. No. 59001-00, is used to update the Odyssey with the latest Hach chemistry methods. This module can be used to update any number of instruments.
2. The Advanced Software Update Module, Cat. No. 59123-00, is used to update the Odyssey with advanced software features, such as scanning, time-course and multiple wavelength analysis. The Advanced Software Update Module is single-use: it can only be used to update one instrument.

Updating the Odyssey

Use the following procedure to update the Odyssey with either module.

1. Turn on the instrument.
2. From the main menu, select **System Checks** then select **Instrument Update**.
3. The instrument will prompt the user to plug the update module into the instrument's serial port. Plug in the module and touch **OK**.
4. The instrument will load the contents of the update module in about two minutes.
5. When prompted, remove the update module from the serial port and touch **OK**. Touch **Return**.

Error Messages

Obsolete Module Error

This error occurs with the Hach Program Update Module (Cat. No. 59001-00) when the chemistry methods contained in the update module are older than those already in the instrument. No update is made when this error occurs.

Module Error

This error indicates that the Hach Advanced Software Update Module (Cat. No. 59123-00) has already been used to upgrade an Odyssey. No additional upgrades can be made with this module.

Components

Description	Cat. No.
Hach Programs Update Module	59001-00
OR	
Advanced Software Update Module.....	59123-00
Instruction Sheet, Update Module.....	59130-00



FOR TECHNICAL ASSISTANCE, PRICE INFORMATION AND ORDERING:

In the U.S.A. – Call toll-free 800-227-4224
 Outside the U.S.A. – Contact the HACH office or distributor serving you.
 On the Worldwide Web – www.hach.com; E-mail – techhelp@hach.com

HACH COMPANY
 WORLD HEADQUARTERS
 Telephone: (970) 669-3050
 FAX: (970) 669-2932