**Sonda 50 60**

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
<tbl\_r cells="2" ix="3" maxcspan="1" maxrspan="1" used

1

50 60


1

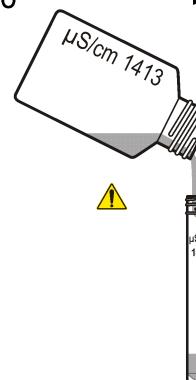2


3

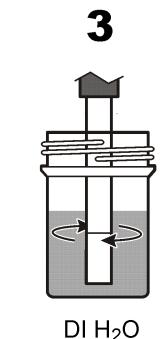4


5

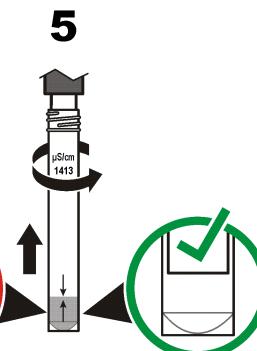


2

4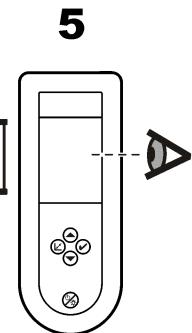

3

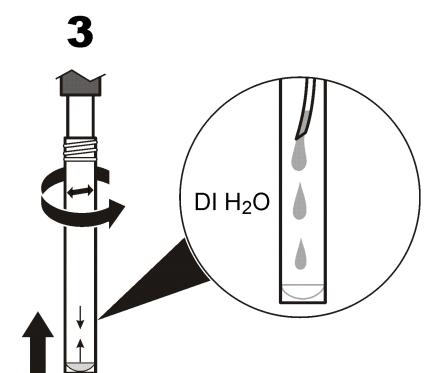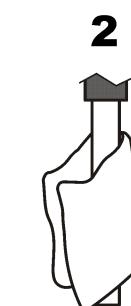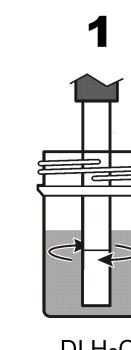EN

**50 60 probe**

Refer to the meter manual for more information.

**Caution**

**⚠ Personal injury hazard. Glass components can break. Handle with care to prevent cuts.**

**⚠ Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.**

**Introduction**

This probe is a 2-pole platinum cell with a built-in temperature sensor for conductivity measurements in general aqueous samples.

**Technical information**

Cell constant	Approximate 1 cm <sup>-1</sup>
Range	0.2 µS/cm to 200 mS/cm
Temperature	Pt 1000
Electrode	Platinum (3x)
Minimum depth	25 mm (1 in.)
Operating temp	0–80 °C (32–176 °F)
Limits	Dirty samples, samples not compatible with polycarbonate or glass.
Connector type	MP-5

**Preparation**

First time use: put the probe in ethanol for 15 s. Rinse with deionized water.

**Connect the probe****Caution**

**⚠ Do not turn or twist the probe connector.**

**1 Calibration****Caution**

**⚠ Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.**

**⚠ Rinse sensor thoroughly between standards.**

For conductivity, use 1-point calibration and select the standard nearest to the expected measurement range.

**2 Measurement**

Continuous: push ✓ 2 times.

Change the parameter: push and hold ✓ during stabilization.

**3 Maintenance and storage**

Do not rub the electrode surface or use an abrasive cleaner.

Cleaning: soak the probe in a solution that dissolves the debris: Inorganic debris: dilute HCl or chromic acid. Organic debris: solvent that is compatible with the probe body. Rinse with deionized water.

**Troubleshooting**

Be sure that the correct probe type is used. Do not hold the tubes by hand to prevent temperature changes.

**Warranty**

1 year for manufacturing faults only. Damage from use is not covered.

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